

# How to install Labview 2020

Note: Following steps are common for Windows 32 bit as well as 64 bit.

1. Download the Unzip Folder from [here](#).
2. Extract the Unzip Folder.
3. Open the unzipped folder and double click on Install.exe to start the installation process.
4. Accept the License agreement.
5. Click on Next to Install NI Package Manager.
6. Select the additional items which you would like to install.
7. Once Again Accept the License agreement.
8. Pop up window comes for NI Update Service. Click on Yes, if you would like the service to check for the updates periodically.
9. For Activation, please contact computer center to Shri Sagar/ Shri Vinit on 022 22782071/2264 respectively.
10. Note your Computer ID. You can open the Labview Application.  
Under Help--->Activate Labview Components--->Launch.
11. NI License Manager will launch. Click on Computer Information. It will display Computer ID.